



# DNA & Biological Evidence Collection

The analysis of DNA is similar to fingerprints in trying to determine a match. With identifying features, DNA can either link or eliminate a suspect from a crime as well as link crime scenes together. The most frequent forms of DNA found at the crime scene are: blood, semen, hair, saliva, skin cells, mucous, perspiration, urine, fingernails, etc. The more we know how to use and collect DNA, the more powerful a tool it becomes. This section contains tools for proper DNA and biological evidence collection.

Evidence	Possible location of DNA on evidence	Source of DNA
Baseball Bat or similar weapon	Handle, end	Sweat, Skin, Blood, Tissue
Hat, Bandana, or Mask	Inside	Sweat, Hair, Dandruff
Eyeglasses	Nose or Ear Pieces, Lens	Sweat, Skin
Facial Tissue, Cotton Swab	Surface Area	Mucous, Blood, Sweat, Semen, Ear Wax
Dirty Laundry	Surface Area	Blood, Sweat, Semen
Toothpick	Tips	Saliva
Used Cigarette	Cigarette Butt	Saliva
Stamp or Envelope	Licked Area	Saliva
Tape or Ligature	Inside/Outside Surfaces	Skin, Sweat
Bottle, Can, or Glass	Sides, Mouthpiece	Saliva, Sweat
Used Condom	Inside/Outside Surfaces	Semen, Vaginal or Rectal cells
Blanket, Pillow, Sheet	Surface Area	Sweat, Hair, Semen, Urine, Saliva
"Through-and-through" Bullet	Outside Surface	Blood, Tissue
Bite Mark	Person's Skin or Clothing	Saliva
Fingernail, Partial Fingernail	Scrapings	Blood, Sweat, Tissue

From the brochure "What Every Law Enforcement Officer Should Know About DNA Evidence", developed under an award from the Office of Justice Program's National Institute of Justice.

## Custom Sexual Assault and Biological Evidence Collection Kits



As a manufacturer, we have full capabilities to create custom kits for the proper collection of DNA and other biological evidence. We work with jurisdictions and organizations around the world to produce kits built to their exact specifications and provide expertise and guidance on best practices to ensure the highest quality up-to-date products.

Contact us today to discuss your needs and receive a tailored end to end solution delivered with world-class customer service.



Often Bought Together:  
Indicates items often purchased in conjunction with other items.

## Sexual Assault Kits

### Sexual Assault Victim Evidence Collection Kit

Recently updated with input from SAFE nurses, this kit is for adult and adolescent examinations and contains all the necessary forms and medical tools to conduct a thorough exam.



#### VEC200A Contents:

- 1- Kit Box, 7" x 10 1/2" x 2 1/4", printed in blue

1- Set Forensic Medical Examination Forms (12 pgs. Total)

1- Step 1a Foreign Material Bag

2- Drape Sheets

2- Step 1B Clothing Bags

1- Step 1C Underwear Bag

1- Step 1D Tampon/Sanitary Napkin Bag

1- Step 2 Debris Collection Envelope containing paper bundle

1- Step 3 Oral Swabs Envelope containing swabs, and swab boxes

1- Step 4 Buccal Swabs Envelope containing swabs, and swab box

1- Step 5 Bite Mark & Breast Swabs Envelope containing swabs, and swab box

3- Step 6 Additional Swabs Envelope containing swabs, and swab box

1- Step 7 Left Hand DNA Collection Envelope containing polyester tipped applicators, swab box, and paper bundle

1- Step 8 Right Hand DNA collection envelope containing polyester tipped applicators, swab box, and paper bundle

1- Step 9 Pubic Hair Combing Envelope containing, paper towel, and black comb
- 1- Step 10 Mons Pubis Swabs Envelope containing swabs, and swab box

1- Step 11 Thigh/External Genitalia Swabs Envelope containing swabs, and swab box

1- Step 12 Perineal Swabs Envelope containing swabs, and swab box

1- Step 13 Vaginal/Cervical Swabs Envelope containing swabs, and swab box

1- Step 14 Penile/Scrotum Swabs Envelope containing swabs and swab box

1- Step 15 Anal/Perianal Swabs envelope containing swabs, and swab box

1- Step 16 Rectal Swabs Envelope containing swabs, and swab box

1- Step 17 Known Blood Sample Envelope containing filter paper

1- Zip Lock Pouch

1- Kiss Marker, Item

1- Kiss Marker, Ruler Item

22- Police Evidence Seals

1- Distilled Water, 1 oz.

1- Swab Box Labels

1- Biohazard Label

1- FDA Sheet

No. VEC200A

Sexual Assault Victim Kit

### Domestic Violence/Strangulation Evidence Collection Kit

Working with Sexual Assault Forensic Examiners (SAFE) across the nation, we developed a Domestic Violence and/or Non-Fatal Strangulation kit. It includes the necessary Forensic/Medical report forms, anatomical drawings, photography log documents, evidence collection envelopes, and measuring tools; all in one place for your convenience. We received feedback from SAFE nurses throughout our development process to ensure you have one of the best kits on the market for your needs.



#### SDV100 Contents:

- 1- Kit Envelope, 9" x 12"

1- Forms Envelope containing:

1- Set Domestic Violence Reporting Officer's Form, 3 pgs.

1- General Information Consent Form, 1 pg.

1- Strangulation History & Event Questions Form, 2 pgs.

1- Domestic Violence Strangulation Exam Forensic Medial Report Forms, 11 pgs.

1- Photography List, 1 pg.

1- Clothing Bag

1- Underwear Bag

1- Debris Collection Envelope containing paper bundle
- 1- Fingernail Swabbings – Right Hand Envelope containing swabs, swab box, and paper bundle

1- Fingernail Swabbings – Left Hand Envelope, containing swabs, swab box, and paper bundle

3- Dried Secretions Envelopes containing swabs, and swab boxes

1- Control Swab Envelope containing swabs, swab box

1- Evidence Seal

1- Measuring Tape

1- Wound Mapping Markers

1- Distilled Water, 1 oz.

1- Biohazard Label

1- FDA Sheet

No. SDV100

Domestic Violence/Strangulation Kit



**Pricing Information:**  
For current pricing visit  
www.sirchie.com, or  
call 800.356.7311

**Often Bought Together:**  
Indicates items often  
purchased in conjunction  
with other items.

Southwestern Sexual Assault  
Evidence Collection Kit

The Texas 100A kit has been designed to meet the specifications of the newly implemented TX Sexual Assault Evidence Tracking System. Created for Texas State Agency use only, the TX100A has all of the components necessary to track the progress of the kit, as well as the status for the victim. The Kit includes a tracking barcode for the state and for the survivors of sexual assault to view the history and current status of the evidence collected by entering the tracking number into a special portal.

No. TX100A Southwestern Sexual Assault Kit

Drug Facilitated Sexual Assault Kits

These kits are used for the collection of blood and/or urine samples from a victim of sexual assault who may have ingested a substance, possibly causing an impairment or incapacitation related to that sexual assault.



DFSABU100 Kit Contents:


- 1- Kit Box, 5.75" x 4.75" x 2.5" with chain of custody label affixed to top, and sealed via integrity seal
- 1- Instruction Sheet/FDA Insert on blue stock, 5.5" x 8.5"
- 1- Information Form, 8.5" x 11", 2-part NCR paper
- 1- 100ml Urine Collection Bottle with PS Liner
- 1- Bubble Pack Bag (double), containing 2 each 10 ml Grey Top Blood Collection Tubes
- 1- Ziplock Bag, 6" x 8", containing Liquid Absorbing Sheet
- 1- Multi Sample Needle, 20g x 1.5
- 1- Tube and Needle Holder with Safety Device
- 1- PVP Prep Pad
- 3- Specimen Security Seals
- 1- Biohazard Label
- 1- Kit Box Shipping Seal

No. DFSABU100 Drug Facilitated Sexual Assault Kit (Blood & Urine)

No. DFSAU100 Drug Facilitated Sexual Assault Kit (Urine)

Pediatric Sexual Assault Kit

Developed with input from SAFE nurses and the state of Alaska, this sexual assault kit was developed to address the unique needs presented in the examination of a possible sexual assault victim who has not yet entered puberty.

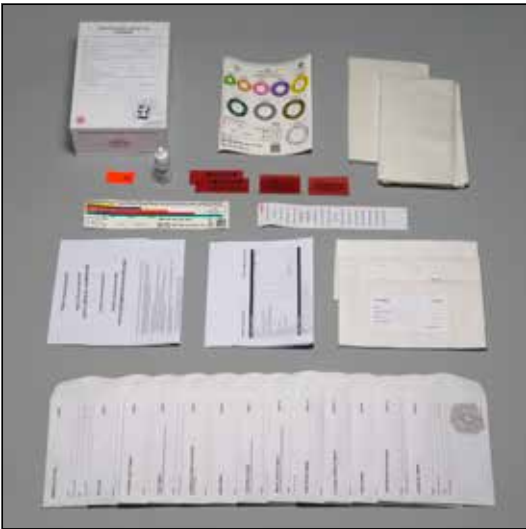

PEDSA100 Contents:

- 1- Kit Box, 7" x 10 1/2" x 3"
- 1- Consent Form
- 1- Photo Documentation Guidelines Form
- 1- Instruction Form
- 1- Step 1 Forensic History Forms
- 1- Step 2 Medical History Forms
- 1- Step 3 Evidence Collection Log Form
- 1- Step 4A Foreign Material Bag
- 2- Step 4B Clothing Bags
- 2- Step 4C Underwear/Diaper Bags
- 2- Drape Sheets
- 1- Step 5 Debris Collection Envelope containing paper bindles
- 1- Step 6 Head Hair Envelope containing paper bindle and black comb
- 1- Step 7 External Mouth Swabs Envelope containing swabs, and swab box
- 1- Step 8 Oral Swabs Envelope containing swabs, and swab box
- 1- Step 9 Known DNA Sample Envelope containing swab, and swab box
- 1- Step 10 Hand Swabs Envelope containing swabs, and swab box
- 1- Step 11 Fingernail Swabs Envelope containing swabs, and swab boxes
- 1- Step 12 Miscellaneous Swabs Envelope containing swabs, and swab boxes
- 1- Step 13 Pubic Mound Swabs Envelope containing swabs, and swab box
- 1- Step 14 External Genitalia Swabs envelope containing swabs, and swab box
- 1- Step15 Penile Swabs Envelope containing swabs, and swab box
- 1- Step16 Perianal/Buttocks Envelope containing swabs, and swab box
- 1- Step 17 Anorectal Swabs Envelope containing swabs, and swab box
- 1- Kiss Marker, item
- 1- Kiss Marker, rings item
- 17- Integrity Seals
- 2- Police Evidence Seals
- 1- Distilled Water, 1 oz.
- 1- Swab Box Labels
- 1- Biohazard label

No. PEDSA100 Pediatric Sexual Assault Kit

Suspect Sexual Assault Kit

Establish culpability or innocence by linking the perpetrator and the victim through DNA with this kit. All necessary components, with clear labeling and step-by-step organization help simplify the collection procedure.



SEC200 Contents:

- 1- Kit Box, 7" x 10 1/2" x 2 1/4" printed in red
- 1- Photo Documentation Guidelines form
- 1- Instructions/FDA form
- 1- Step 1 Release Form
- 1- Step 2 Patient History Forms
- 1- Step 3A Foreign Material Bag
- 2- Step 3B Clothing Bag
- 2- Step 3C Underwear Bag
- 2- Drape Sheets
- 1- Step 4 Debris Collection Env w/paper bindle
- 1- Step 5 Dried Secretions Env w/swabs & swab boxes
- 1- Step 6 Head Hair Combing Env w/paper towel & black comb
- 1- Step 7 External Mouth Swabs Env w/swabs & swab box
- 1- Step 8 Oral Swabs Env w/ swabs & swab box
- 1- Step 9 Known DNA Samples Env w/ swabs & swab box
- 1- Step 10 Fingernail Swabs Env w/swabs & swab box
- 1- Step 11 Hand Swabs Env w/ swabs & swab box
- 1- Step 12 Pubic Hair Combing Env w/paper towel & comb
- 1- Step 13 Penile Swabs Env w/swabs & swab box
- 1- Step 14 Scrotal Swabs Env w/swabs & swab box
- 1- Step 15 Anal/Perianal Swabs Env w/ swabs & swab box
- 1- Step 16 Rectal Swabs Env w/swabs & swab box
- 1- Evidence Collection Log Form
- 1- Kiss Marker, Item
- 1- Kiss Marker, Rings Item
- 1- Swab Box Labels
- 1- Distilled Water, 1 oz
- 1- Police Evidence Seals, 20/pk
- 1- Biohazard Label

No. SEC200 Suspect Sexual Assault Kit

Suspect Evidence  
Collection Kit (Buccal  
Swab/Finger Stick)

Collect evidentiary samples from a sexual assault suspect. Instructions guide through collecting and preserving evidence from clothing, hair samples, buccal swabs & blood.



SEC100BF Kit Contents:

- 1- Kit Envelope, 7.5" x 10.5"
- 1- Instruction Sheet/FDA Form
- 1- Step 1 Underpants Bag
- 2- Step 1 Outer Clothing Bag
- 1- Step 2 Foreign Material Collection Env w/ paper bindles, scrapers, swabs & swab box
- 1- Step 3 Pubic Hair Combing Env w/ paper towel & comb
- 1- Step 4 Pulled Pubic Hairs Env
- 1- Step 5 Pulled Head Hairs Env
- 1- Step 6 Buccal Swab Env w/ swabs & swab box
- 1- Step 7 Known Blood Sample Env w/ 21G Haemolance Lance, alcohol swab, gauze pad, adhesive bandage & bloodstain card
- 1- Step 8 Anatomical Drawing Sheet
- 1- Evidence Seal
- 1- Biohazard Label

No. SEC100BF Suspect Collection Kit (Buccal Swab/Finger Stick)

Portable Swab  
Drying Kit

This Portable Swab Drying Kit allows you to dry up to 36 swabs at a time. Simply collect sample swabs; place them in the included Swab Drying Racks; place racks into the kit case, and close the lid. Once the swabs are dry, place them into the included swab boxes for storage, and transport.



SDK100 Contents:

- 1- Plastic Carrying Case
- 5- Swab Drying Racks
- 10- Swab Boxes
- 10- Cotton-Tipped Applicator Swab Packs (2/pack)

No. SDK100 Portable Swab Drying Kit

Wound Map Marker Kit

These markers are a simple but highly effective method to establish the size, placement, orientation, and identification of wounds. Using this kit will help you document and track wounds and help keep your photos organized. Marker sheets are also available in individual packs: circular wounds, map markers, rulers, stick figures, and ellipse wound guide.



No. WMK80 Wound Map Marker Kit, 8 Markers, 10 ea.

No. WMK80CR Circular Wound Guides, 10 ea.

No. WMK80ER Ellipse Wound Guides, 10 ea.

No. WMK80CR9 Wound Guide Circle Scales, 1-9, 10 ea.

No. WMK80R25 Wound Guide Rulers, 10 ea.

No. WMK80SF Wound Guide Stick Figures, 10 ea.

No. WMK80MM Wound Guide Map Markers, 1-14, 10 ea.

No. WMK80M50 Wound Guide Map Markers, 1-50, 10 ea.

No. WMK80M100 Wound Guide Map Markers, 51-100, 10 ea.

**Pricing Information:**  
For current pricing visit  
www.sirchie.com, or  
call 800.356.7311

**Often Bought Together:**  
Indicates items often  
purchased in conjunction  
with other items.

Forensic Swab Dryer

The **New and Improved** Sirchie Forensic Swab Dryer is manufactured with a **stainless steel cabinet** that is designed to stand up to frequent cleanings between drying sessions, especially when harsh chemicals like chlorine bleach solution is used. It's compact, dries swabs efficiently and can be secured to protect chain of custody while swabs are drying.

Use in forensic laboratory or forensic work area or anywhere where evidence swabs are taken, providing a safe, controlled environment to dry swabs for processing later. It holds up to 6 swab-drying racks (up to 72 swabs in total) and will dry them to greater than 99% dry in about an hour.

The Forensic Swab Dryer uses HEPA filtration with axial fans pulling air through HEPA filters with additional filtering on the exhaust to eliminate biological contaminants from entering the room. It has a digital timer with two ranges: 1-999 seconds or 1-999 minutes. An alarm/chime will sound to alert you when the set cycle is complete.

A latch is included with non-tamper tags that can be used to secure the door to protect chain of custody while drying.

FSD100 Features:

- Compact size: suitable for table/benchtop use and lightweight for easy movement as needed
- Low energy consumption
- Low fan noise
- Modern appearance with interior cool white lighting
- Easy to set drying time with two ranges: 1-999 seconds or 1-999 minutes
- Digital times that emit audible (2 beeps) and visual indicator (displays END) when selected drying time is complete

Forensic Swab Dryer includes:

- 1- Drying cabinet
- 2- HEPA intake-side filters
- 1- Fabric exhaust filter

1- Starter kit with:

- 2- swab boxes that hold 12 swabs each
- 5- blue security ties
- 5- red security ties

No. FSD100	Forensic Swab Dryer, 110V AC
No. FSD100220	Forensic Swab Dryer, 220V AC
No. FSD01	Replacement Exhaust Filter, 1 ea.
No. FSD02	Replacement HEPA Filters, 2 ea.



Swabbing Kits

Buccal Swab Kit

With the Buccal Swab Kit, non-medical personnel can collect a buccal cell sample for DNA testing by swabbing the inside of the subject's cheek. Swabs are then placed inside the provided envelope and forwarded to the laboratory for analysis.



BSK100 Contents:

- 1- BSK1000 Kraft Mailing Envelope 7.5" x 10.5" (19cm x 26.7cm)
- 1- BSK101 Kit Shipping Seal
- 1- BSK1001 White Kit Env. 6" x 9" (15.2cm x 22.9cm) with Factory Seal
- 1- BSK1002 Instruction Sheet
- 1- BSK103 Information Card with Attached Swab Envelope
- 1- KCP124 Pouch, Silica Gel
- 1- KCP125 Zip-Top Bag 5" x 8' x .004" (12.7cm x 20.3cm x 4 mil)
- 1- KCP160 Sterile, Cotton-Tipped Swabs, 2 pk.
- 2- KCP164 Evidence Seals, Red, "Warning! Police Seal, Do Not Remove"

No. BSK100	Buccal Swab Kit
------------	-----------------

Uni-Swab Kit

Avoid cross-contamination of swabs after collecting samples by using swabs with applicator shields. Swabs are individually sterilized.



US100 Contents:

- 1- US001B Kit Box; Dimensions: 9.25" x 7.25" x 1.375"
- 20- US001 Uni-Swabs consisting of:
  - 1- KCP160 Applicator
  - 1- AS24 Swab Shield
- 1- TX101 Swab Box
- 2- ECT2A Biohazard Seals
- 1- US04 Chevron Pouch
- 1- US02 Instruction Sheet

No. US100	Uni-Swab Kit
-----------	--------------

Child DNA Identification Kit

This kit has been designed to collect information, fingerprints, photographs, and a DNA specimen from a child. In the event the child goes missing, this kit is presented to law enforcement to assist in finding and identifying the child.

DNACID100 Contents:

- 1- Kit Envelope, 6" x 9", sealed with tamper tape
- 1- Information Sheet/FDA, 11" x 14"
- 1- Saliva Swabs Envelope, 3 ½" x 6 ½"
- 1- Pack of Cotton Tipped Swabs, 2 each
- 1- Security Seal
- 1- Pre-inked Fingerprint Sheet
- 1- Photographs (CD/DVD), Glassine Envelope, 5 ½" x 5 ½"

No. DNACID100	Child DNA Identification Kit
---------------	------------------------------

Fingernail Scrapings, Swabbings & Clippings Kit

This kit contains the items necessary for the collection of fingernail scrapings, swabbings and clippings for both hands as well as an evidence seal to ensure evidence integrity.

FSSC100 Contents:

- 1- 6" x 9" Kit Envelope, sealed with tamper tape
- 1- Instruction Sheet/FDA Insert
- 1- Left Hand Scraping Envelope w/paper bindle & plastic scraper
- 1- Left Hand Clippings Envelope w/paper bindle
- 1- Right Hand Scraping Envelope w/paper bindle & plastic scraper
- 1- Right Hand Clippings Envelope w/paper bindle
- 1- Swab Box labeled "Right Hand Finger Swabs"
- 1- Swab Box labeled "Left Hand Finger Swabs"
- 1- Package, Cotton-Tipped Applicator Swabs (2/pkg)
- 1- Distilled Water, 3cc
- 1- Evidence Seal
- 1- Fingernail clipper

No. FSSC100	Fingernail Scrappings, Swabbings & Clippings Kit
-------------	--



Blood & Urine Collection Kits

Blood Specimen Collection Kit

Designed for medical personnel, this kit has everything to collect a whole blood sample for DNA testing, preserve it, and send to the laboratory.

BSC50 Contents:

- 1- White Corrugated Mailer, 5" x 4" x 2.5"
- 1- Instruction Sheet/FDA Insert
- 1- Information Card
- 1- Iodine Prep Pad, Medium
- 1- Multi-Sample Needle, 2G x 1.25" w/Luer Adapter
- 1- Blood Collection Tube, 7ml with EDTA
- 1- Specimen Security Seal
- 1- Absorption Sheet, 2.5" x 4"
- 1- Zip-Top Evidence Bag, 4" x 6" x .004"
- 1- Biohazard Label, 1" x 1.25"
- 1- Kit Shipping Seal
- 1- Gauze Pads, 12-Ply, 2" x 2"
- 1- Adhesive Backed Bandage
- 1- Bubble Pouch, 4" x 5.5"
- 1- Ink Strip, 1" x 5"
- 1- Needle, 21 ga. x 1.25" w/Luer Adapter

No. BSC50	Blood Specimen Collection Kit
-----------	-------------------------------

Blood/Urine Specimen Collection Kit

This kit contains all the necessary components for taking blood samples and urine collection for substance abuse testing. Follow the enclosed step-by-step instructions and fill out all documentation. Package the specimens for analysis using the enclosed plastic bag with absorbent material and protective pouches. Seal inside the crush resistant white container and deliver to the laboratory.

BUK200 Contents:

- 1- White Corrugated Mailer, 5" x 4" x 2.5"
- 1- Instruction Sheet/FDA Insert
- 1- Consent Form
- 1- Police Officer's Report
- 1- Blood/Urine Collection Report
- 1- Iodine Prep Pad, Medium
- 1- Multi-Sample Needle, 2G x 1.25" w/Luer adapter
- 2- Blood Collection Tube, 10ml
- 1- Needle, 21 ga. x 1.25" w/Luer adapter
- 3- Specimen Security Seal
- 1- Absorption Sheet, 2.5" x 4"
- 1- Zip-Top Bag, 6" x 8" x .002"
- 1- Evidence Seal
- 1- Biohazard Label, 1" x 1.25"
- 1- Bubble Pouch, 4" x 5.5"
- 1- Urine Specimen Jar, 100ml

No. BUK200	Blood/Urine Specimen Collection Kit
------------	-------------------------------------





**Pricing Information:**  
For current pricing visit  
www.sirchie.com, or  
call 800.356.7311

**Often Bought Together:**  
Indicates items often  
purchased in conjunction  
with other items.

Blood Card  
Collection Kit

This kit contains all the necessary components to collect 100 dried blood samples for the long term preservation of DNA for up to 50 years (based on US DOD standards), and is a great low-cost alternative to preserving dried blood samples with expensive, traditional DNA collection cards.

BBC100E Contents:

- 1- White Corrugated Mailer, 5 3/4" x 4 1/4" x 3 1/2", integrity sealed
- 100- Manila Coin Envelopes, 3 1/2" x 2 1/4"
- 100- Subject/Case Identifier Labels for envelopes, 1 1/8" x 2"
- 100- Blank Blood Cards, 2 1/4" x 3"

No. BBC100E Blood Card Collection Kit



Blood Alcohol  
Collection Kit

Our comprehensive blood alcohol collection kit is designed to provide the safe and efficient collection, handling, and mailing of blood specimens to the lab.

BSC100A Contents:

- 1- FDA Manufacturer's Insert
- 1- Police Officer's Report
- 1- Instruction Sheet
- 1- Consent Form
- 1- Blood Collection Report
- 1- Cardboard Tube Holder/Protector
- 1- Needle, 21 ga. x 1.25" with Luer Adapter
- 2- Blood Collection Tubes, 10ml with Gray Top
- 1- Iodine Prep Pad, Medium
- 1- Liquids Absorption Sheet
- 1- Multi-Sample needle: 2G x 1.25" w/Luer Adapter
- 1- Tourniquet, Latex
- 1- Zip-Top Bag, 6" x 9" x 0.004
- 1- Evidence Seal
- 2- Blood Tube Integrity Seals
- 1- Biohazard Label
- 1- FDA Manufacturer's Insert
- 1- Cardboard Kit Mailer

No. BSC100A Blood Alcohol Collection Kit



Blood Alcohol/Urine  
Specimen Collection  
Kit

This kit is ideal for departments with multi-sample requirements. It offers the same safety and efficiency as our Blood Alcohol Collection Kit but with the added collection tools for a urine specimen.

BUK100A Contents:

- 1- Instruction Sheet
- 1- Police Officer's Report (Blood)
- 1- Police Officer's Report (Urine)
- 1- Consent Form for Blood & Urine
- 1- Collection Report for Blood & Urine
- 1- Cardboard Blood Tube & Needle Protector
- 2- Blood Collection Tubes, 10ml with Gray Top
- 1- Needle, 21 ga. x 1.25" (3.2cm) w/Luer adapter
- 1- Tourniquet, Latex
- 1- Needle Holder
- 1- Urine Specimen Container, Sterile, 120ml
- 1- Iodine Prep Pad, Medium
- 2- Liquids Absorption sheet
- 2- Zip-Top Bag, 6" x 9" x 0.004"

No. BUK100A Blood Alcohol/Urine Specimen Collection Kit



Postmortem Blood  
Urine Specimen Kit

The Post-mortem Blood Urine Specimen kit is designed specifically for gathering specimens from a deceased body. Everything you need in one place to take blood samples or collect urine for testing.

PMBU100 Contents:


- 1- Kit box with biohazard label
- 1- FDA product insert
- 1- Blood tube and urine bottle holder with security flap
- 1- 100ml Urine specimen bottle
- 2- 10ml blood collection tubes
- 1- 16 gauge and 4" needle
- 1- Ziplock bag with a liquid-absorbing sheet
- 3- Tamper-Evident Specimen security/identification seals
- 1- Tamper-evident kit box shipping seal for resealing outer box after specimen collection

No. PMBU100 Postmortem Blood Urine Specimen Kit



Trace Evidence Collection Kit

Designed to collect dry trace evidence, use these tools to preserve trace evidence at a crime scene.

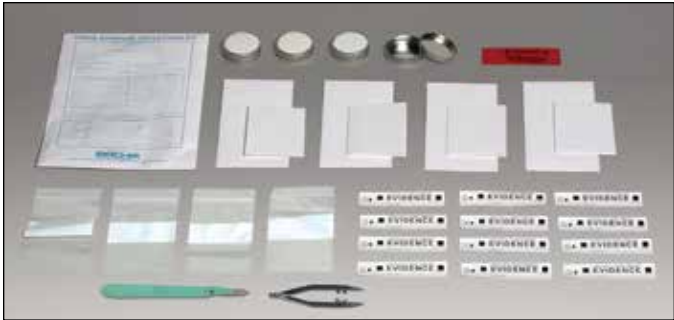

TEC100 Contents:

- 1- Factory-Sealed White Kit envelope, 7.5" x 10.5"
- 4- Evidence Collection Containers, 1 oz
- 4- Write-Block Zip-Top Evidence Bags, 3" x 5"
- 4- White Evidence Collection Envelopes, 3.5" x 6.5"
- 4- Unfolded Paper Bindle, 8.5" x 11"
- 1- Sterile Scalpel
- 12- Evidence Seals for Evidence Containers, 3.25" x 1"
- 1- Red Evidence Seal

No. TEC100 Trace Evidence Collection Kit

DNA Collection Kits

DNA Specimen Collection Belt Kit

Keep supplies close at hand with this kit that clips directly on your belt. It contains everything you'll need for DNA Specimen Collection.



DNA15 Contents:

- 16- KCP307 Unprinted Swab Boxes
- 8- KCP208 Sterile Cotton Swabs, 2 pack
- 16- KCP206 Black on Orange, Biohazard Labels, 1" x 1.25"
- 4- KCP247 Sterile Water, 3ml
- 1- Molded Plastic, Belt-Style Case, 2.25" x 4.25" x 4"

No. DNA15 DNA Specimen Collection Belt Kit

Crime Scene DNA Evidence Collection Kit

You should always be prepared to collect potential sources of DNA and this kit has everything you need in one place. The most frequent forms of this valuable evidence are: fresh blood, semen, hair roots, or clothing stained by physiological fluid.



DNA100 Contents:

- 1- Kit Box, 10.25" x 6.75" x 2.25"
- 2- Evidence Identification Strips
- 1- FDA Manufacturer's Insert
- 1- Information Sheet
- 5- Evidence Collection Envelopes
- 1- Undergarment Bag
- 1- Debris Collection Bag
- 2- Outer Clothing Collection Bags
- 1- Tape Strip, Red, Vinyl, 19"
- 1- Distilled Water, 1 oz.
- 4- Evidence Bag Integrity Seals
- 2- Blood Collection Tubes, 7ml w/EDTA
- 1- Syringe w/Needle, Disposable, 10cc
- 1- Pair of Tweezers, Plastic, 4.75"
- 5- Sterile, Cotton-Tipped Swabs, 2pk
- 2- Evidence Seals
- 1- Liquids Absorption sheet
- 1- Mailer, Styrofoam, Two-Tube
- 5- Swab Box, Ventilated
- 4- Biohazard Label
- 5- Gauze Pads, 12-Ply, 2" x 2"
- 1- Zip-Top Evidence Bag, 6" x 9"
- 6- Latex Powder-Free Gloves

No. DNA100 Crime Scene DNA Evidence Collection Kit

**Pricing Information:**  
For current pricing visit  
www.sirchie.com, or  
call 800.356.7311

**Often Bought Together:**  
Indicates items often  
purchased in conjunction  
with other items.

DNA/Biological Evidence Collection Kit

Stored in a rugged case with a sliding drawer, you'll easily find everything you need to collect, store, and mark DNA and other biological evidence at crime scenes.



DNA200 Contents:

- 10- Sterile Cotton-tipped Applicator Swabs, 2/pk
- 20- Swab Box, printed
  - 1- Swab Drying Rack
  - 2- Clear Polystyrene Jars w/lid, 1 oz.
  - 2- Clear Polystyrene Jars w/lid, 2 oz.
  - 2- Clear Polystyrene Jars w/lid, 4 oz.
  - 1- LED Penlight w/2 ea. AAA batteries
- 20- White Coin Env, printed, 3.5" x 6.5"
- 5- Mask, disposable tie
  - 1- Straight Hemostat, 5.5" stainless steel
  - 1- Curved Scissors, 4.5"
  - 1- Spiral Bound Notebook, 4" x 6", 50 sheets
- 20- "Biohazard Handle with Gloves" Seals, 1" x 4"
- 10- Integrity Seals
  - 1- Felt Tip Marker, extra fine, black
- 4- Scalpel, round blade, disposable #10, sterile
- 10- Evidence ID Labels, 2.75" x 1.75"
  - 1- Distilled Water, 1 oz.
- 3- Black Nitrile Gloves, Pair
- 5- Preprinted Evidence Bags, 7" x 4.5" x 13.75"
  - 1- Brochure "What Every Law Enforcement Officer Should Know About DNA Evidence"
  - 1- High-Impact Copolymer Carrying Case w/handle

No. DNA200 DNA/Biological Evidence Collection Kit

Applicator Swabs & Accessories

We offer a variety of swabs and accessories to meet your needs for collecting evidence at crime scenes:

Sterile Cotton Applicator Swabs and Swab Boxes

Ideal for the collection of trace liquid evidence, such as blood and other bodily fluids. Includes swab boxes so you can mark important information on the box.



No. KCP160C	Sterile Cotton Swabs, 6" (2 pk), 100 ea.
No. KCP160M	Sterile Cotton Swabs, 6" (2 pk), 1000 ea.
No. KCP208C	Sterile Cotton Swabs, 3" (2 pk), 100 ea.
No. CC1094C	Preprinted Swab Boxes, 6", 100 ea.
No. CC1094M	Preprinted Swab Boxes 6", 1000 ea.
No. KCP196C	Plain Swab Boxes, 6", 100 ea.
No. KCP196M	Plain Swab Boxes, 6", 1000 ea.
No. KCP307C	Plain Swab Boxes, 3", 100 ea.

Fab-Swab DNA Collection Applicator

This convenient, reclosable, dry transport device is designed for reliable sample and evidence collection. The patented tube has a breathable filter label that covers the swab to allow drying while protecting against contamination.



No. FSWAB50 Fab-Swab DNA Collection Applicator, 50 ea.

Cap-Shure® Sterile Forensic Evidence Collection Swabs

These swabs are a unique crafted method for collecting small quantities of very fragile evidence. Whether it's blood, saliva, latent print residue or other physiological fluids, these Cap-Shure® Swabs ensure the integrity of the samples.



No. CP100	Cap-Shure® Sterile Swabs, 100 ea.
No. CP500	Cap-Shure® Sterile Swabs, 500 ea.
No. CP50PL	Cap-Shure® Sterile Swabs, 50 ea.
No. KCP254C	Vented Swab Boxes, 3/4" x 3/4" x 6 1/4", 100 ea.

Applicator Shields

To prevent swab cross-contamination after collecting samples, as well as during shipment to the laboratory, use these sturdy applicator shields.



No. AS24 Applicator Shields, 24 ea.

Sterile Round Foam-Tipped Applicators

These over-sized, foamed-tipped swabs have a large surface area that permit the collection of larger samples than common, cotton-tipped swabs.



No. FTS50W Sterile Round Foam-Tipped Applicators, 50 ea.

Sterile Omni-Swabs

The Sterile Omni-Swab's unique design uses an absorbent multi-layered paper collection and a non-reactive material that allows DNA extraction for analysis. Additionally, the swab's head is ejectable and fits into an extraction tube after collection, making it easily transported to the lab.



No. SOS100 Sterile Omni-Swabs, 100 ea.

Cuticle Stick

Manufactured in a controlled environment and individually wrapped, these double-beveled, contaminant-free cuticle sticks are ideal for collecting evidence.



No. CSTICK125 Cuticle Stick, 125 ea.

HydraFlock® Sterile Swabs

HydraFlock swabs produce superior sample collection, rapid elution, and increased test sensitivity by combining a unique microstructure with innovative fibers.



No. HFS50 HydraFlock Sterile Swabs, 50 ea.

Purflock® Sterile Swabs

Made with nylon fibers, these swabs have the ability to better absorb organisms at the collection site and during transport.



No. PFS50 Purflock Sterile Swabs, 50 ea.

CollectEject™ Swabs


These ejectable-head collection swabs make collecting and processing DNA samples efficient and easy. Collected samples are stabilized & preserved with solid desiccation with improved DNA yield, and the ejectable swab head helps prevent cross contamination.



No. XXXXX CollectEject™ Swabs, XX ea.

GenSwab™ Collection Swabs

GenSwab™ is designed for noninvasive collection, preservation, shipment, and long-term storage of DNA samples at room temperature. Collected samples are stabilized & preserved with solid desiccation with improved DNA yield, and can be stored at ambient temperature for 24 years without degradation. Each swab is individually barcoded for sample management.

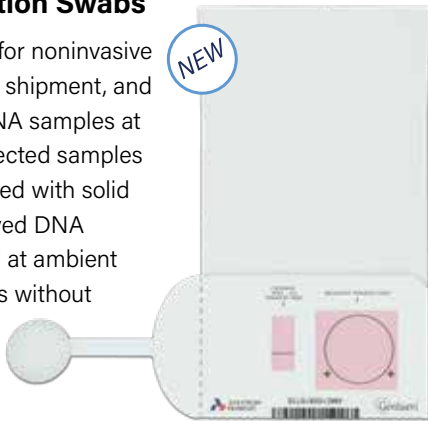

No. XXXXX GenSwab™ Collection Swabs, XX ea.

Swab Drying Rack

Use the Swab Drying Rack to reduce the risk of cross contamination while applicators are drying. It is constructed from 20 mil (0.020") thick board material with a built-in inner board shelf that holds up to 12 applicators upright and straight. The rack top is preprinted with collection location descriptions. This Swab Drying Rack can also be used to air dry swabs from any type of specimen collection.



No. SDR100C Swab Drying Rack, 100 ea.



**Pricing Information:**  
For current pricing visit  
www.sirchie.com, or  
call 800.356.7311

**Often Bought Together:**  
Indicates items often  
purchased in conjunction  
with other items.

EZ-DRY™ Swab boxes and Evidence/Integrity seals

The EZ-DRY™ Swab Box is one of the best ways to collect and package cotton swabs used to collect biological evidence—they dry at room temperature. Once the collection swabs have been packaged in the swab box, use our EZ-DRY™ Swab Box Evidence/Integrity Seals. Fill out the necessary information on the label and then apply it to the swab box, overlapping the ends in order to form the integrity seal.



No. KCP248C	EZ-Dry Swab Boxes, 100 ea.
No. SB102C	EZ-Dry Swab Box Evidence/Integrity Seals, 100 ea.

EZ-Dry™ Swab Box Kit

This EZ-Dry™ Swab Box Kit is the best way to collect, package, label and seal cotton swabs without the possibility of sample switching or secondary contamination, while providing a way to maintain a continuous chain of possession. Room temperature is sufficient to dry swabs with EZ-DRY™— no freezer or swab dryer is needed.



- SB1000 Contents:**
- 1- SB1001 Kit Envelope
  - 5- KCP248 EZ-DRY Swab Boxes
  - 5- KCP160 Sterile Swabs, Pack of 2
  - 1- Distilled Water, 3ml
  - 5- SB102 Evidence/Integrity Seals
  - 1- Instruction Sheet
  - 5- KCP206 Biohazard Labels, 1" x 1 ¼", Black on Orange

No. SB1000	EZ-Dry Swab Box Kit
------------	---------------------

Sterile Cotton Swabs

Easy-to-use cotton tip swab with a long handle to collect evidence. Available in 2-packs of 100 or 1000 with wooden handles or with plastic handles in a 2-pack of 100.



No. PUR8062WC	Sterile Cotton Swabs, wood handle, 2pk, 100 ea.
No. PUR8062WM	Sterile Cotton Swabs, wood handle, 2pk, 1000 ea.
No. PUR8062PC	Sterile Cotton Swabs, plastic handle, 2pk, 100 ea.

DNA-Free Specimen Collection

DNA Collection Cards

These collection cards are treated with a proprietary chemistry intended to prevent environmentally induced degradation. They allow long term preservation of DNA from biological fluids stored at ambient temperature. Use *non-indicating* cards for blood sampling, and *indicating* cards (that turn color from pink to white) for saliva, buccal cells and urine.



No. DNAG1	DNA Card, 1-Spot (non-indicating), 100 ea.
No. DNAG2	DNA Card, 2-Spot (non-indicating), 100 ea.
No. DNAG4	DNA Card, 4-Spot (non-indicating), 100 ea.
No. DNAG1C	DNA Card, 1-Spot (indicating), 100 ea.
No. DNAG2C	DNA Card, 2-Spot (indicating), 100 ea.
No. DNAG4C	DNA Card, 4-Spot (indicating), 100 ea.

DNA-Free Swabs

Certified to be DNA-Free, these applicators are made with high-quality USP grade cotton.



No. FDNA100	DNA-Free Swabs, 100 ea.
-------------	-------------------------

DNA-Free Ultra Flocked Swabs

DNA-Free swab with a plastic handle and standard tip.



No. FDNA350	DNA-Free Ultra Flocked Swabs, 50 ea.
-------------	--------------------------------------

SecurSwab™ 2

These DNA-Free swabs come with barcodes to provide the end-user with control in labeling the sample and maintaining chain-of-custody control.

No. SWB100N	SecurSwab 2, 100 ea.
No. SWB100B	SecurSwab 2 w/barcode, 100 ea.

SecurSwab™ Duo-V

This is a complete system that includes components for collection, preservation & drying, protection, tracking & transport. Available with or without barcodes.

No. SDV100N	SecurSwab Duo-V, 100 ea.
No. SDV100B	SecurSwab Duo-V w/barcode, 100 ea.

Presumptive Seminal Fluid Identification

The Acid Phosphatase Test

In sexual assault cases, you can save some time with a presumptive non-destructive chemical test to identify possible seminal stains.

The Acid Phosphatase Test is one of those presumptive tests and it's considered to be one of the most conclusive, non-spermatozoa tests for the presence of semen. The reagents for this test are sealed in glass ampoules to ensure integrity and long life. You should conduct a preliminary exam with longwave UV lights or an alternate light source, such as BLUEMAXX (BM300)

The reagents should not be used directly on the suspect stain. Using the contact method, a filter paper wetted with distilled water is pressed against the stain, and the paper is then tested. This preserves the stain for later confirmatory and possible DNA testing.

Seminal ID AP Test

The Acid Phosphatase (AP) Test is used for the presumptive identification of semen and seminal fluid. Sirchie now offers this field test in an easy-to-use, rugged pouch format. Each box includes 10 tests, filter paper, and distilled water.



Easy-to-Use:

1. Simply dampen a piece of the included filter paper with a few drops of the provided distilled water.
2. Press the filter paper to the suspected stain.
3. Place the filter paper in the pouch, and close the pouch with the safety clip.
4. Gently squeeze the ampoule in the pouch and agitate.
5. A color change to a purple is a positive result for the presence of semen or seminal fluid.

No. PSID100	Seminal ID AP Test, 10 tests
-------------	------------------------------

**Pricing Information:**  
For current pricing visit  
www.sirchie.com, or  
call 800.356.7311

**Often Bought Together:**  
Indicates items often  
purchased in conjunction  
with other items.

Master Seminal Fluid Test Kit for Detection of Acid Phosphatase

This is a complete Master Field and Lab Kit that enables the investigator to perform contact testing and verification of suspected seminal stains. Components of the various reagents are sealed in glass vials to eliminate contamination and assure unlimited storage life. Applicators are disposable. Because of the specificity of the test, only minute quantities are required for identification and stains up to four years in age have responded positively. The contact transfer technique is used to verify the presence of semen eliminating contamination probability.



- SF298 Contents:**
- 8- 288DP Pipettes, Disposable with Capillary and 3ml Bulb
  - 2- 288FP Contact Filter Paper, 1" x 2"
  - 2- DISW2 Bottles of Distilled Water, 2 oz
  - 10- EIL011 Evidence Identification Labels, "Crime Evidence"
  - 3- KCP172 Glass Vials with Cap, 3dr
  - 15- KCP217 Cotton Balls
  - 8- SF298S1 S1 Reagent
  - 8- SF298S2 S2 Reagent
  - 8- SF298S3 S3 Reagent
  - 1- SF298C Black, Molded Copolymer Carrying Case with Folding Handle & Custom Insert, 14" x 10.75" x 5.75"

No. SF298 Master Seminal Fluid Test Kit

RSID Semen Detection Field Test Kits

This presumptive test for semen is great to use in the field as a confirmatory test for human semen. The test is specific for human semenogelin antigen. There is no cross reactivity with other body fluids or animal semen. Strip test results are complete in 10 minutes. Each kit includes 10 test units and protocol.



Each individual RSID™ Semen test includes:

- 1- Test kit (reagent, pipette, swab, test cassette)
- 1- Zip-Top bag
- 1- Sterile water, 3ml
- 1- Documentation envelope

No. IDSE200 RSID Semen Detection Field Test Kit, 10 tests

RSID Saliva Detection Field Test Kit

This presumptive test for saliva is great to use in the field as a confirmatory test for human saliva. The test is specific for human salivary (amyA) antigen. There is no cross reactivity with blood, semen, urine, vaginal secretions, or menstrual blood. Strip test results are complete in 10 minutes. Each kit includes 10 test units and protocol.



Each individual RSID™ Saliva test includes:

- 1- Test kit (reagent, pipette, swab, test cassette)
- 1- Zip-Top bag
- 1- Sterile water, 3ml
- 1- Documentation envelope

No. IDSA300 RSID Saliva Detection Field Test Kit, 10 tests

RSID Bodily Fluid Detection Field test Kit

This kit combines 4 presumptive blood tests, 4 presumptive semen tests, and 2 Presumptive saliva tests and protocols.



No. IDBF400 RSID Bodily Fluid Detection Field Test Kit, 10 tests

Specimen Collection Tubes and Accessory Products

Sterile & Distilled Water

Pre-moisten applicator swabs when collecting dry specimens with sterile or distilled water. Avoid over-saturating swabs.



**Often Bought Together:**  
(KCP160C) Sterile Cotton Applicator Swabs

No. KCP247C Sterile Water, 3ml vial, 100 ea.

No. DNA1009 Non-Sterile Distilled Water, 1 oz. bottle, 1 ea.

Blood Collection Tubes

**Gray stopper blood collection tubes:** 10 ml, 16 x 100mm with no interior coating contains 20mg potassium oxalate and 100mg sodium fluoride. **Lavendar stopper blood collection tubes:** 7 ml, 13 x 100mm contains K3 EDTA 15% solution, 0.081 ml, 12.15 mg. **Red stopper blood collection tubes:** 10 ml, 16 x 100mm, silicone coated.



No. KCP176C Gray Stoppered Blood Tube, 10 ml, 100 ea.

No. KCP122C Lavendar Stoppered Blood Tube, 10 ml, 100 ea.

No. CC110C Red Stoppered Blood Tube, 10 ml, 100 ea.

Padded Blood Tube Containers

These 1" x 4 1/2" containers are large enough to hold one 10ml blood tube. Protector pad included.



**Often Bought Together:**  
(KCP176C) Gray Stoppered Blood Tube, (KCP122C) Lavendar Stoppered Blood Tube, (CC110C) Red Stoppered Blood Tube

No. BTC100C Blood Tube Container and Plug Insert, 100 ea.

Blood Tube Containers and Caps

These containers are large enough to hold one 10ml tube providing extra protection.



**Often Bought Together:**  
(KCP176C) Gray Stoppered Blood Tube, (KCP122C) Lavendar Stoppered Blood Tube, (CC110C) Red Stoppered Blood Tube

No. BTC200C Blood Tube Containers & Caps, 100 ea.

Bubble Pouch Tube Protectors

These bubble pouch tube protectors have a peel and seal flap that keeps the tubes secure in the pouch.



**Often Bought Together:**  
(KCP176C) Gray Stoppered Blood Tube, (KCP122C) Lavendar Stoppered Blood Tube, (CC110C) Red Stoppered Blood Tube

No. KCP253C Bubble Pouch Tube Protectors, 100 ea.

Liquid Absorption Sheets

Ideal anywhere where bodily fluids are being collected, these sheets are also perfect to ship along with specimen containers during transport.



**Often Bought Together:**  
(KCP176C) Gray Stoppered Blood Tube, (KCP122C) Lavendar Stoppered Blood Tube, (CC110C) Red Stoppered Blood Tube

No. KCP113C Liquid Absorption Sheets, 2 1/2" x 4", 100 ea.

Tube Protector Pads

For added protection, wrap blood collection tubes in tube protector pads before inserting them in containers or bubble pouches.



**Often Bought Together:**  
(KCP176C) Gray Stoppered Blood Tube, (KCP122C) Lavendar Stoppered Blood Tube, (CC110C) Red Stoppered Blood Tube

No. KCP305C Tube Protector Pads, 100 ea.

Toluidine Blue O Solution

This clinically proven staining agent, Toluidine Blue O, is used during exams of sexual assault victims to reveal and visualize any lesions caused during an assault that are not visible without an enhancement agent. Unlike single-use popules products, our dye has 15 applications per bottle, enabling multiple locations to be tested for lesions without stopping to open single-test packaging for each test area. It speeds up the exam for the victim, which reduces their stress. Sold in a 1 oz. (30ml) pre-mixed bottle.



No. TBD100 Toluidine Blue O Solution